# **Cole-Parmer**®

## Cole-Parmer® Paraffin Wax Dispenser - WD-200

- 'Wax on tap' for histology and pathology labs
- Over-temperature cut-out at 120°C
- Non-drip heated tap to prevent blockages
- Anti-microbial coating (Ag+) to inhibit growth of bacteria
- Slimline design to maximise bench space







### **Technical Specification**

Specification	WD-200		
Capacity	7.5 L		
Material	Stainless Steel		
Temperature	Ambient to 70°C		
Dimensions (d x w x h)	500 x 181 x 455mm		
Weight	7.0 kg		

## Cole-Parmer<sup>®</sup> Paraffin Wax Dispenser - WD-200

The Cole-Parmer<sup>®</sup> paraffin wax dispenser is designed for on-demand delivery of molten wax. It heats wax to a liquid state and provides convenient 'wax on tap' for Histology and Pathology laboratories. The paraffin wax dispenser has a large 7.5 L capacity tank, for melting up to 6 kg of pelletised wax at a time.

The innovative design of the instrument ensures that it takes up minimal bench space in the laboratory. The wax dispenser has digital temperature control which provides accurate temperature control of the molten wax, and is coupled with an ultra-fast heating system for rapid melting of pelletised wax.

An excellent feature is that the wax dispenser tap is heated via a dedicated heating system which prevents blockages and ensures an even flow of wax ondemand. Delivery of molten wax is via a non-drip lever tap with three settings (off, on-demand pull dispense and continuous flow). The inner tank has a filter screen (0.5 mm mesh) fitted to prevent coarse particles from blocking the delivery tap.

The wax dispenser is controlled via an intuitive digital control panel. The set temperature can be controlled using the up and down arrows on the control panel. The heater indicator will illuminate to show heater activity. The wax dispenser is designed to melt bulk quantities of pelletised wax in as quick a time as possible. When set at  $65^{\circ}$ C the instrument will take ~90 minutes to fully melt 5 kg of pelletised wax, with significant quantities of molten wax available for use after as little as 30 minutes.

The equipment is protected by a miniature circuit breaker and a temperature cut out. If an over-temperature condition occurs, the wax dispenser will automatically turn off at 120°C.

### **Ordering Information**

Description	Ordering Number	Series No.	Model No.	Legacy Sku.
Wax Dispenser 230V	85230-07	WD-200	WD-200D	MH8524
Wax Dispenser 120V	85230-08	WD-200	WD-200D-120	MH8524X1

## coleparmer.com

USA: +1.800.323.4340 +1.847.549.7600 Canada: +1.800.363.5900 China: +86.21.5109.9909 France: +33 (0) 1486 37800 Germany: +49 (0) 9377 92030 India: +1.800.266.1244 Italy: +39 (0) 1313 89513 UK: +44 (0) 1480 272279 All other countries: +1.847.549.7600 Cole-Parmer essentials an antylia scientific company

AZ-EMEA-LT-00245 1121

Find out more at antylia.com